

BL 520 PRO-ACTIVE

according to DIN 13277

Equipment

- 5 aluminium drawers
- Forced-air cooling
- Equipped with castors
- Key switch
- Automatic defrosting
- Condensate evaporation
- Potential-free contact
- Digital temperature display
- Comfortable access
- Antifreeze
- Visual and audible alarm signal, even in the event of power failure
- Minimum/maximum temperature memory

MADE IN GERMANY  0483



Interior may differ visually.

PRO-ACTIVE-Features

- Faster response time in the event of impending faults thanks to proactive alarms
- Data recording can be read out via USB interface using a USB stick and software
- Minimisation of temperature deviations in the cold room
- Maximum storage safety thanks to harmonised components



- **PRO-ACTIVE-Control:** Permanent, proactive monitoring of the performance data and alerts in the case of deviations. World's most accurate temperature control in refrigerated areas thanks to two standard PT-1000 sensors.
- **External housing** made from galvanised sheet steel (rust-proof), with white, anti-scratch powder coating. Length of the power cable: approx. 3 m.
- **Equipped with castors.**
- **Interior made from** smooth aluminium with colourless protective coating. Rails (every 15 mm) for flexible interior arrangement.
- **Interior** can be customised (surcharge may apply). See optional equipment.
- **Extra-thick insulation**, made from high-quality, compression-moulded and environmentally friendly material. Energy saving.
- **Self-closing door** with easy-to-replace plastic magnetic seal frame, lockable.
- **Door stop** on the right-hand side (by default, see illustration), or the left-hand side, can be retrofitted.
- **Forced-air cooling**, with axial blower, switches off automatically when you open the door, ensures a uniform temperature and minimizes temperature deviations.
- **Automatic fast defrosting** thanks to time-limited and temperature-monitored reversal of the refrigerant circuit.
- **Condensate evaporation** in stainless steel heated dish beneath the refrigerator floor.
- **Control panel** protected against tampering by key switch. Membrane keyboard with digital temperature display and minimum/maximum temperature memory.
- **Power failure alarm** (visual and audible alarm), the monitoring unit remains in operation for approx. 30 hours on battery power.
- **Warning functions** with visual and audible alarm signal in the case of temperature deviations or other malfunctions. Door open alarm after 60 seconds.
- **Alarms** can be forwarded via email or SMS using the Kirsch EVOLUTION CLOUD Module (optional).
- **Data documentation** can be read via USB interface with KIRSCH Datanet software.
- **Antifreeze** against sub-zero temperatures.
- **Ventilation-enforced refrigerating machine** hermetically sealed, energy saving, low-noise, easy to service.
- **Low-noise compressor** reduces noise.

Specification

| | |
|--|--------------------------------|
| Capacity | 500 litres |
| Temperature setting | +4 °C |
| Voltage | 220 – 240 V, 50/60 Hz |
| Power consumption | 60 watts |
| Normal consumption | 0.89 kWh/24 h |
| Admissible ambient temperature | from +10 °C to +38 °C |
| Heat emission (max.) | 554 watts |
| Exterior dimensions (including door handle and distance from wall) | w x d x h = 77 x 76 x 195.5 cm |
| Usable dimensions | w x d x h = 56 x 39.5 x 129 cm |
| Exterior dimensions with door open at 90° | w x d = 83 x 144 cm |
| Drawer inner dimensions | w x d x h = 56 x 39.5 x 10 cm |
| Max. load drawer | 50 kg |
| Weight (net/gross) | 144 kg/167 kg |
| Temperature deviation | +/-0.8 °C |
| Noise emission | 47.7 dB(A) |

Optional equipment

KIRSCH refrigerators and freezers offer numerous equipment options that can be retrofitted.



Glass door with lock

Transparency: avoid opening the door unnecessarily to inspect the contents. Lockable.




Additional Shelves

Wire shelves, robust with load capacity up to 40 and 50 kg.



Door coupling

Connects the furniture door to the device.

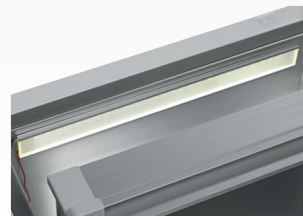

Interior lighting, side wall

Optional, optimal and energy-efficient lighting for the interior using an LED lightstrip.



Decorative frame

Made of stainless steel, allows to cover door with a decorative panel (thickness max. 2 mm).



Condensation tray

For manual draining of condensation water.




Access Port

Simple insertion of an external temperature sensor with integrated cable feed-through.



Levelling feet

Levelling feet instead of castors.

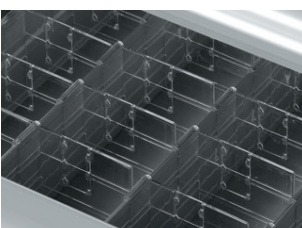

Additional length and cross dividers

For the optimum organisation of your refrigerated goods.



Drawers on runners

Up to eight drawers are possible in total.



Vario Support


Storage shelf on runners.

Accessories for temperature documentation




PC-KIT-NET

Automatic temperature documentation and monitoring via your network (LAN connection).

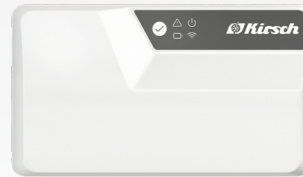

KIRSCH EVOLUTION CLOUD-Module

Maximum storage safety through optimal connectivity and real-time monitoring. *

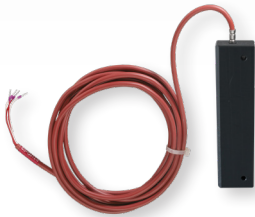

External PT100/PT1000 with cold block

Additional temperature sensor for connection to existing/external monitoring system.



Circular blade temperature recorder

Depending on the model, it is either installed in the machine compartment panel of the unit or integrated into an additional housing attachment.

* For more information about our cloud, please contact us.

Philipp Kirsch GmbH

Im Lossenfeld 14
77731 Willstätt

Telefon: +49 (0) 781 9227-0
Telefax: +49 (0) 781 9227-200
info@kirsch-medical.de

www.kirsch-medical.com



You can find our complete product overview on our website:
www.kirsch-medical.com